

## H&E Protocol Sheet – For Leica Autostainer

(Modified from Leica Protocol)

No.	Station	Reagent	Time min:s	Exact	Dip	Notes
1	Oven	65°C	10	No		
2	1	Xylene	3	No	3	
3	2	Xylene	2	No	3	
4	3	Xylene	2	No	3	
5	4	100% Alcohol	1	No	3	
6	5	100% Alcohol	1	No	3	
7	6	95% Alcohol	1	No	3	
8	Wash 1	H <sub>2</sub> O	1:30	No	3	
9	8	HAEMATOXLIN	2	Yes	3	Change 2000 slides or two weeks
10	Wash 5	H <sub>2</sub> O	1	No	3	
11	9	DEFINE	:35	Yes	3	Change daily
12	Wash 4	H <sub>2</sub> O	1	No	3	
13	10	BLUE BUFFER	:45	Yes	3	Change weekly
14	Wash 3	H <sub>2</sub> O	1	No	3	
15	11	95% Alcohol	1	No	3	
16	12	EOSIN	1	Yes	3	Change 1500 slides or 1- 2 weeks
17	13	100% Alcohol	1	No	3	
18	14	100% Alcohol	1	No	3	
19	16	100% Alcohol	1	No	3	
20	17	Xylene	1	No	3	
21	18	Xylene	1	No	3	
22	22	EXIT				